



User Manual

Alltech customer

What is Alltech® RAPIREAD™?

Alltech RAPIREAD is the latest addition to our integrated system of tools that provide the ultimate protection from the threat of mycotoxins. Suitable for testing individual ingredients and forages, the system highlights real-time mycotoxin contamination on-farm or in the feed mill, providing you with 24/7 online access to detailed results and analysis.

The Alltech RAPIREAD system consists of:

- The Raptor® Integrated Analysis Platform.
- Alltech RAPIREAD Mobile App.
- Alltech RAPIREAD Data Sync Tool.
- Mycotoxin Management Portal.

Accurate results can currently be determined in the field or in the feed mill environment for aflatoxin, ochratoxin, zearalenone, DON, T2/HT-2 Toxins, fumonisin and ergot.

Important: Please read all instructions in this document before proceeding to submit a test request and carry out mycotoxin testing with Alltech RAPIREAD.

Table of Contents

Page 1

- What is Alltech RAPIREAD?

Page 2

- Creating an account on the Mycotoxin Management Portal

Page 4

- Downloading and installing the Data Sync Tool on your PC or laptop

Page 6

- Downloading the Alltech RAPIREAD mobile app on your IOS or Android device

Page 9

- Submitting a mycotoxin test on the Alltech RAPIREAD app

Page 11

- Carrying out the mycotoxin test on the Neogen Raptor device

Page 13

- Uploading test results using the Data Sync Tool

Page 16

- Viewing mycotoxin quick reports on the Mycotoxin Management Portal

Page 18


- Viewing a mycotoxin quick report on the Alltech RAPIREAD app

Page 21

- Overview video and further questions

There are three key steps you should take before submitting your first mycotoxin test request on Alltech RAPIREAD.

Step 1. Create your personalized account on the Alltech Mycotoxin Management Portal

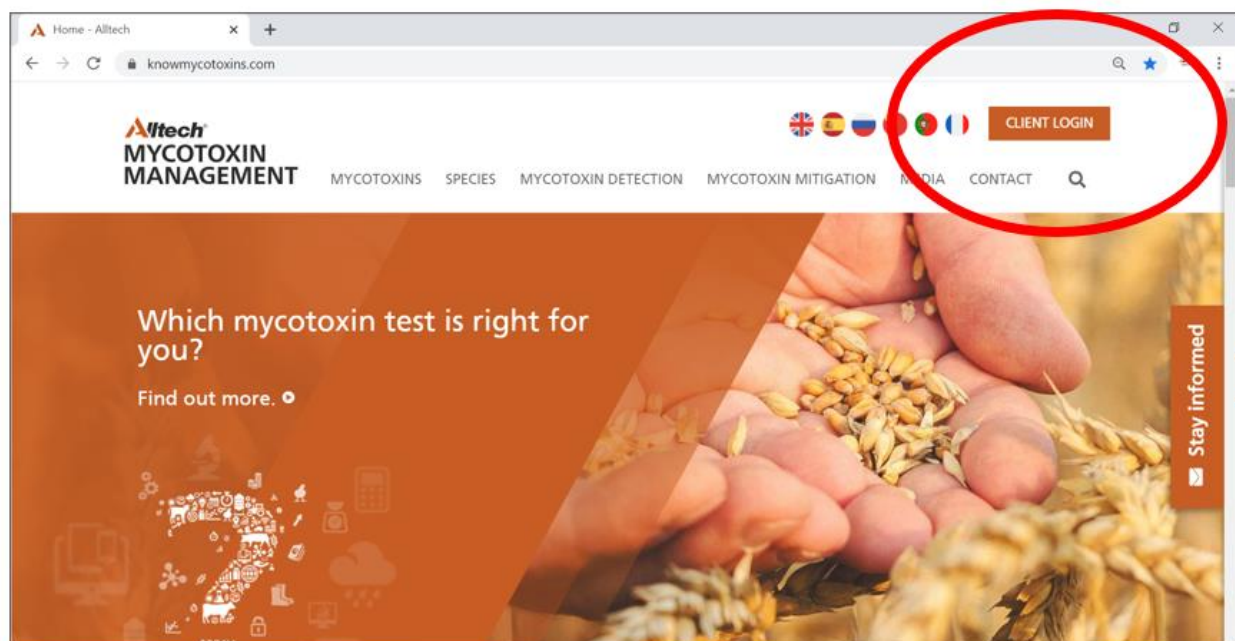

- A. Visit knowmycotoxins.com and click **Client Log In** on the top right of the page
- B. Click on the drop-down in the top right to select country and language
- C. Click on the **Customer Sign Up** on the **User Log In** Panel.
- D. Complete all required fields.


Tip: Contact your office manager before signing up to ensure a quick turnaround of account approval. If you are unsure of who your office manager is please email jteulker@alltech.com.

Note: Please select the level of access you require on the **User Level** drop down tab. **Basic** level is suitable for users who are linked with one account. **Advanced** level is for those who are to be linked to more than one account e.g. nutritionists or quality control managers that have multiple mill associations.


Select **both Alltech 37+ Test and Rapid Test** option on the **Type** drop-down tab.

Step 2. Download and install the Data Sync Tool on your PC or laptop


- A. Log into the Mycotoxin Management Portal at www.knowmycotoxins.com and click CLIENT LOGIN.
- B. Enter your user credentials and Login.
- C. Click the **Data Sync Tool** found on the red header and wait for the setup file named 'publish.zip' to download.

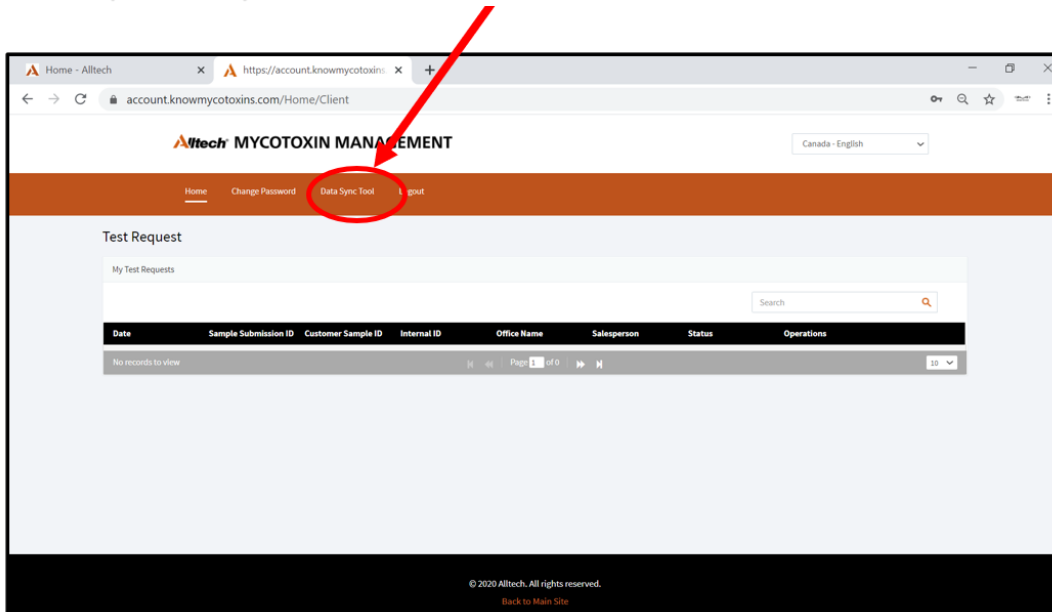

- D. Once downloaded, click the download icon in the top left corner. If a Google Drive warning appears, click **download anyway**.

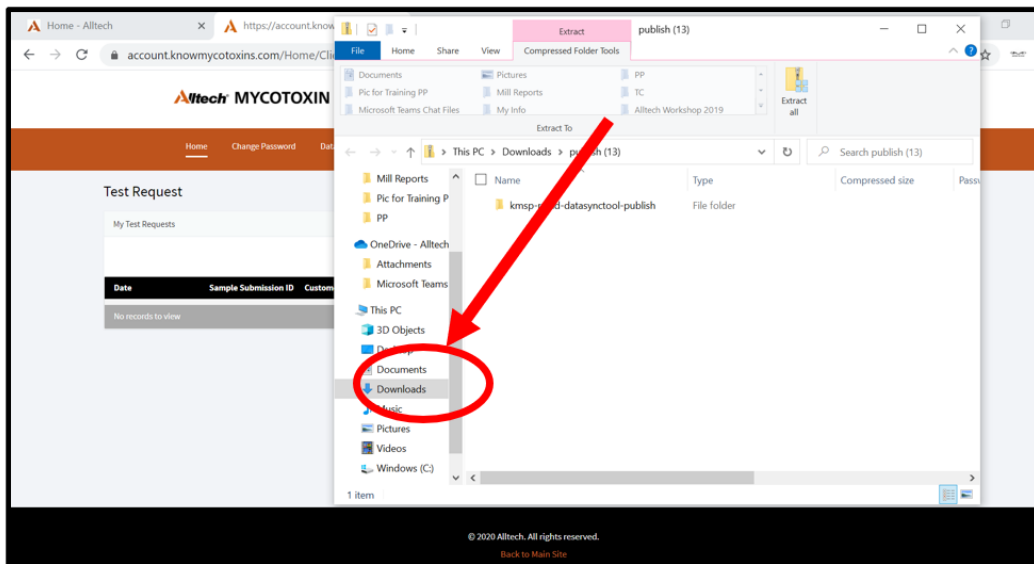

- E. Once the file has downloaded, open the publish folder and double click the setup file. If a windows defender warning appears, click **More Info** and then click **Run Anyway**.

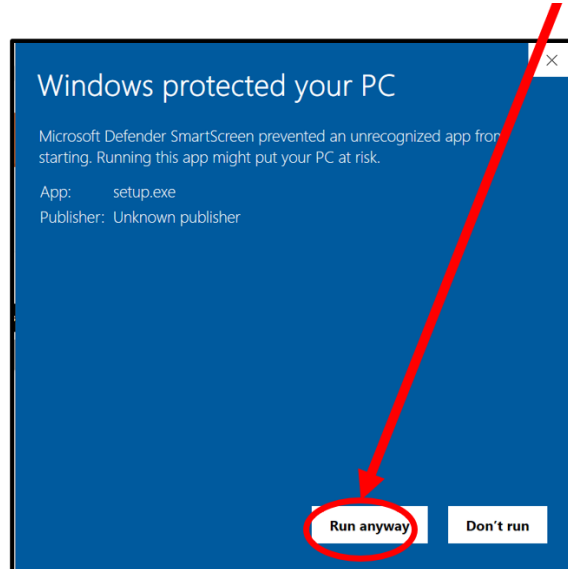


- F. Click **Install** and wait until the program fully installs
- G. The Data Sync Tool should now be installed, and you will be taken to the sign-up screen.
- H. You will now be able to access the Data Sync Tool from your windows desktop under **RAPIREAD Data Sync Tool**.

Signing into the Data Sync Tool:


1. Open the Data Sync Tool on your PC.
2. Enter your username and password. (Same credentials as the Mycotoxin Management Portal)
3. Click Submit.
4. Go to your emails and wait for the verification code email. Copy and paste the verification code into the appropriate fields.
5. Click Sign In.
6. The Data Sync Tool will now open and you can begin to sync files.

Step 3. Downloading the Alltech RAPIREAD Mobile App on your IOS or Android device


- A. In the App or Play Store search **Alltech RAPIREAD**.
 - a. Click GET for IOS App Store and enter your Apple ID and username to install.
 - b. Click Install for Android users in the Play Store to install.

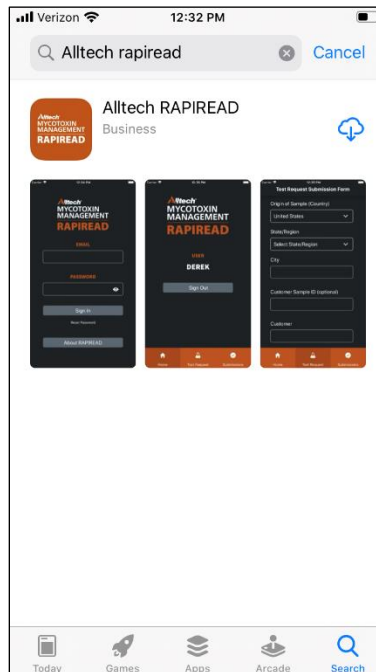


- B. Once installed, open the app and you will be taken to the app's sign in page. Sign in with your portal email and password.

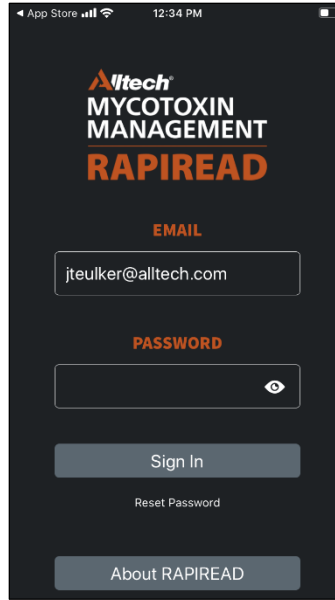

C. A verification code will be sent your email.

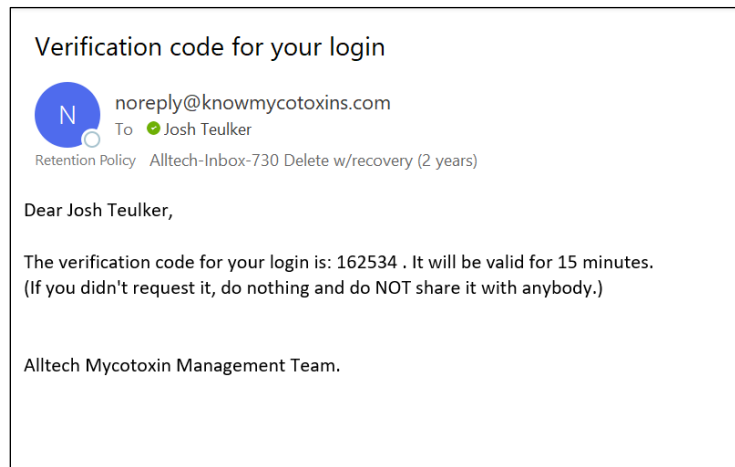

D. Enter the verification code and click **Verify**.

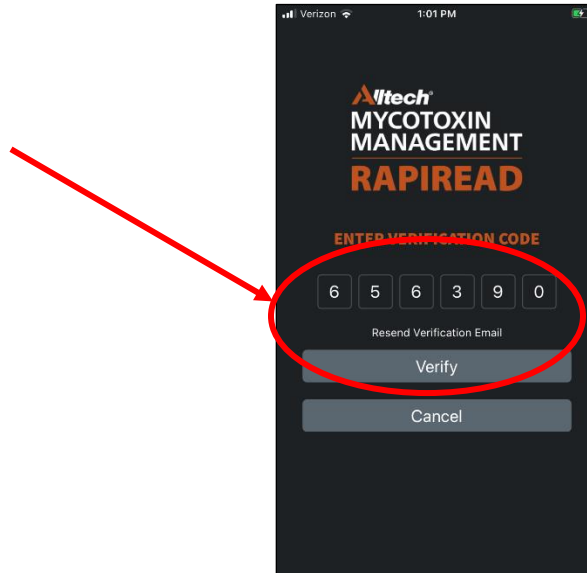

Congratulations! You have now downloaded and installed each of the necessary tools to run mycotoxin tests and view the subsequent results on Alltech RAPIREAD.


Note: You will be required to login every 30 days. The previous login will expire after 30 days and you will require a new verification code when signing in again.

Submitting a mycotoxin test on the Alltech RAPIREAD app

Note: The same Mycotoxin Management Portal login credentials must be used for the Data Sync Tool and the Alltech RAPIREAD App to successfully carry out a test and transfer the information.

Step 1. Create a test request using the Alltech RAPIREAD App

- A. Open the Alltech RAPIREAD App and Login if required.
- B. Click on Test Request tab at the bottom of the screen.

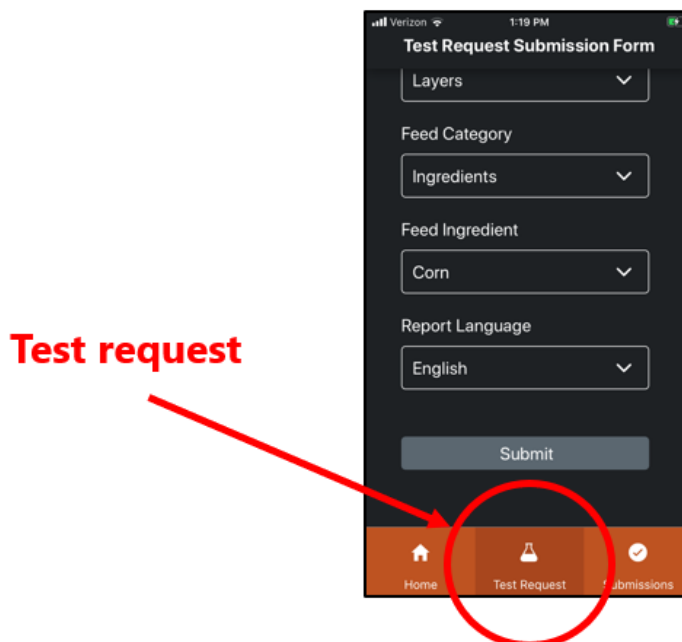

- C. Complete all the required fields. When filling in the 'customer' tab, select the appropriate customer from the list.
 - If you are creating a test for a specific salesperson, select the office the salesperson is associated with. For example: If you are running a test for Nick Adams, you would select the 'Office-UK' account. A 'contact' drop down will appear and then you would select Nick Adams.
 - i. **Tip:** All the fields on the test request tab can be prefilled to save time. For Example: If a test is always going to be ran in a certain State/Region that box can be prefilled.
- D. Click Submit.

- E. Check that all details are correct on the submission page and click **Confirm** to finish.

- F. A QR Code will be generated. This code will be used on the Raptor in future steps.

Tip: If you lose the QR code you created go to the **Submissions** tab on the App and click on the test request. This will bring the QR code back up on the app.


Note: Ensure you have good internet or mobile connection when creating a test request. If connection is interrupted the test request may not work properly.

Step 2. Carrying out the mycotoxin test on the Neogen Raptor device

- A. Follow the guidelines Neogen provide to prepare a sample for testing. These guidelines are found in the booklets provided in each Reveal Quantitative Test Box.



- B. Once the sample is prepared and ready to test insert the appropriate test strip in a cartridge and insert the cartridge into 1 of the 3 ports on the top of the Raptor.
- C. The Raptor will read the test strip and ask to either Enter Sample ID or Scan Sample ID. Choose **Scan Sample ID**.

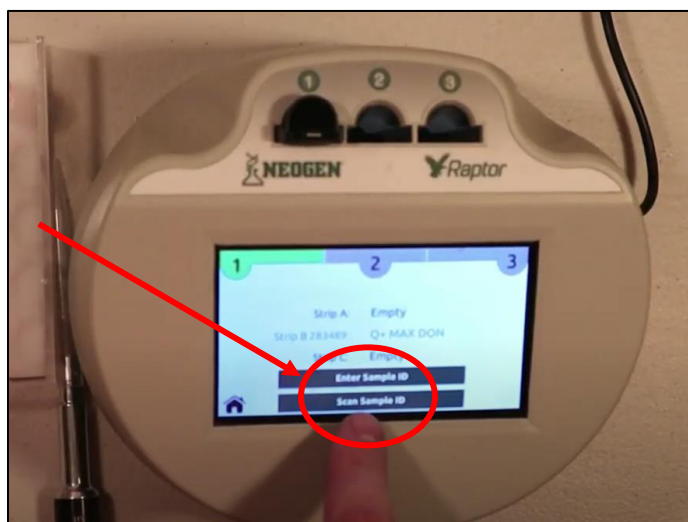

D. Scan the QR Code that was created on the App for the sample being ran.



Tip: Ensure the backlight on your phone is turn to the brightest setting.

E. Continue to follow Neogen's instructions to complete the mycotoxin test.


Note: Each QR Code represents 1 sample. Continue to scan the same QR code if testing for multiple mycotoxins. 3 or more toxin tests are required for all features such as REQ on the quick report.

Step 3. Uploading test results using the Data Sync Tool


- A. Once all mycotoxin tests have been completed for the sample, plug the use the USB cable to connect the Neogen Raptor device to your laptop or PC.
 - a. Note: This step will only work via USB. Network connection will result in an error.




- B. Open and log in to the Alltech RAPIREAD Data Sync Tool if necessary.
- C. On the Raptor, select the computer icon and select **connect with USB** option.

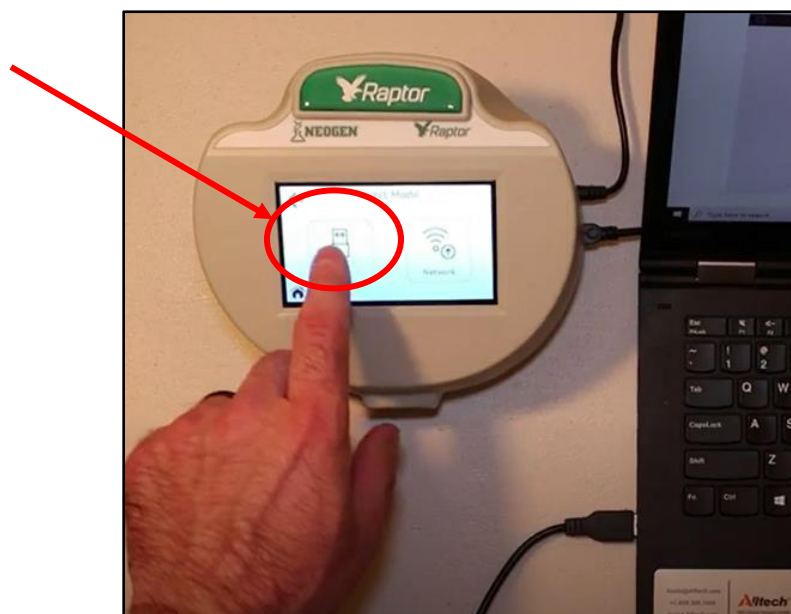

- D. It may take a few minutes for the Raptor to install. Wait until your PC displays the **Device is ready** message before continuing.
- E. Once installed, click the **Connect** button in the right corner of the screen. All files on the Raptor should now display under the **Files** tab.

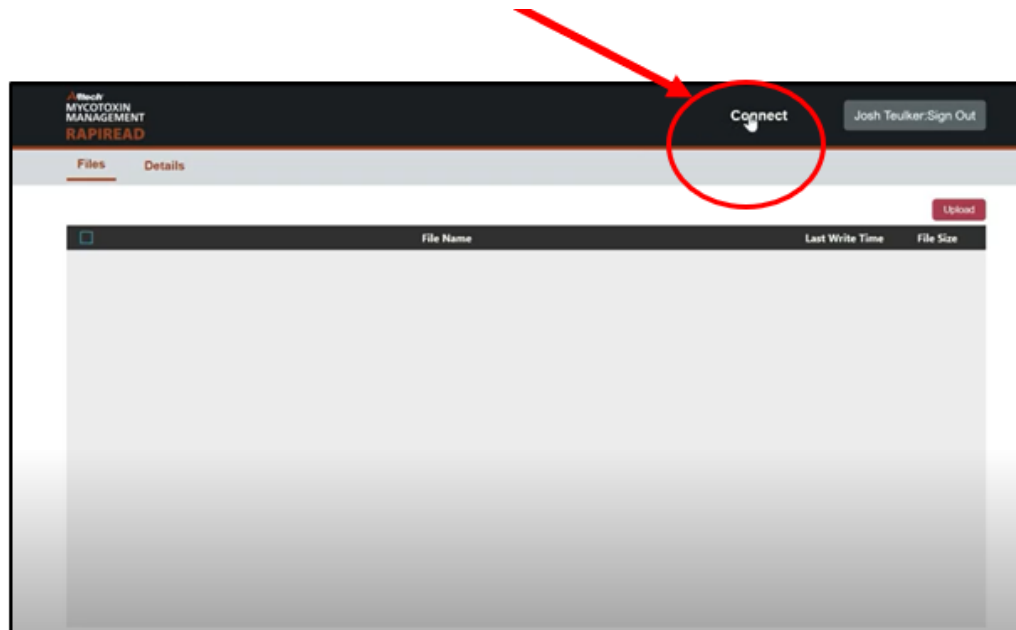


- F. Select the files you want to sync by clicking on the check box on the left side of the file name.

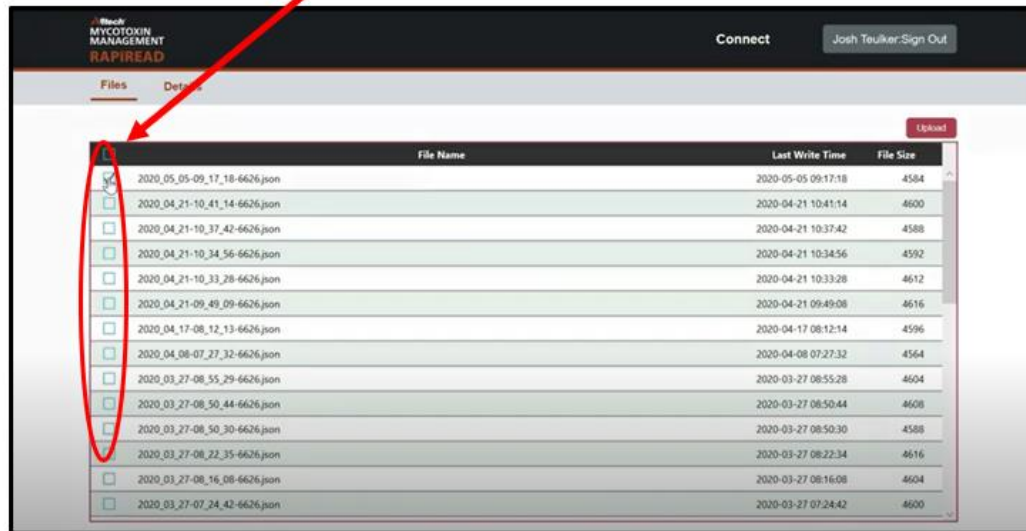

Note: An error will occur if any test is selected that have already been uploaded to Alltech RAPIREAD or does not have a QR Code associated with the sample.


- G. To check the details of a test sample, click the **Details** tab.
- H. Click **Upload** and wait for the files to finish syncing.

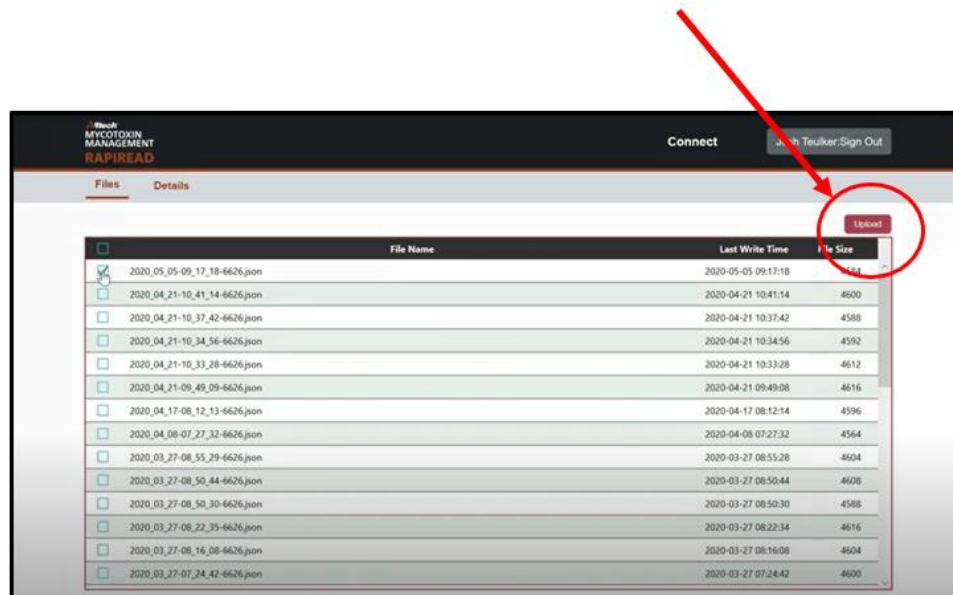


- I. Your report will now be available to view.

Step 4. Viewing mycotoxin quick reports on the Mycotoxin Management Portal


- Visit www.knowmycotoxins.com.
- Click on **Client Log In** and enter your Username and Password.
- Each sample will be located on the home screen.
- Click **View Reports** under the Operations column.

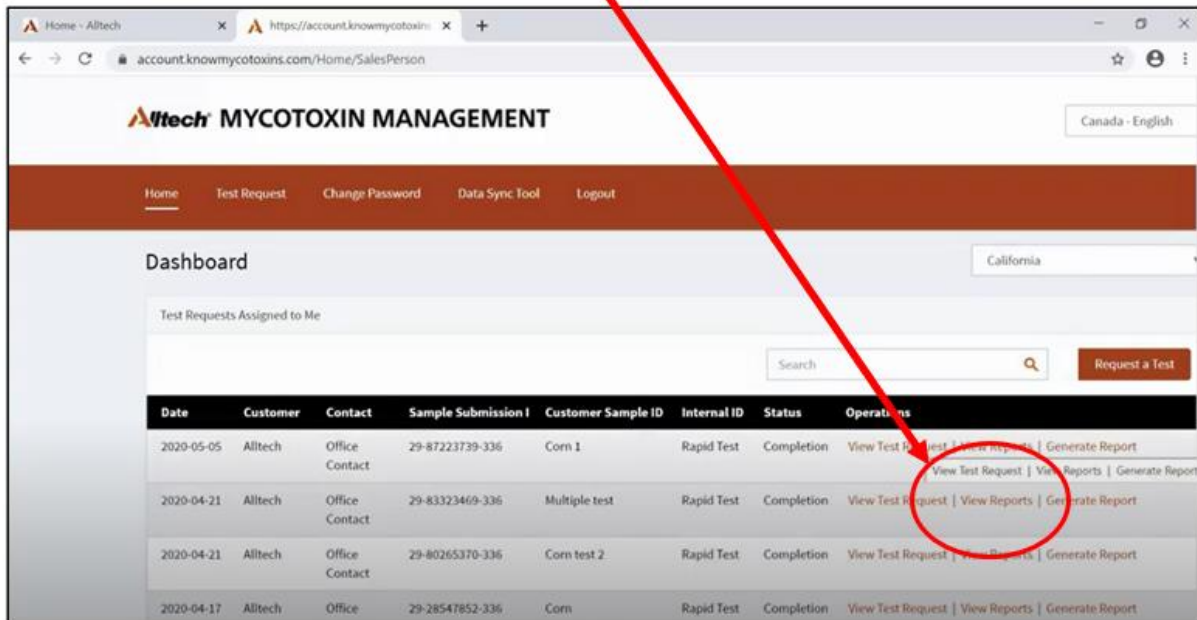

- Scroll to the bottom of the under Final Reports and click **Assessment Report** under the Reports column.

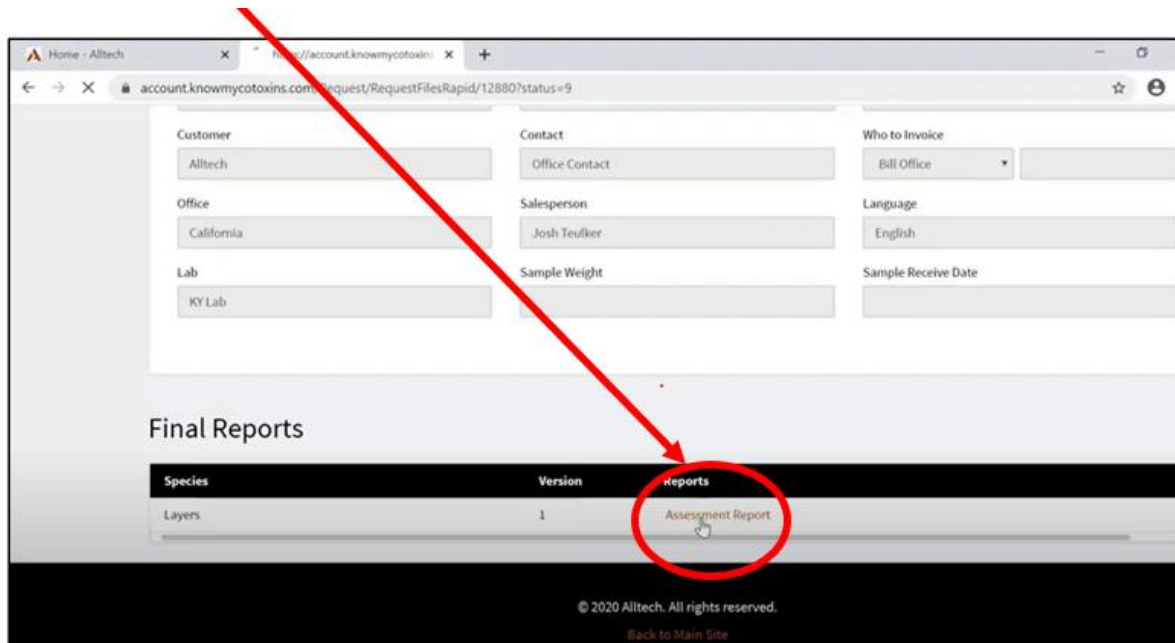


Notes: You can also generate new reports for different species by clicking **Generate Report** under the operations column on the home screen. Then select a new species in the dropdown box and click **Generate Report** in the bottom right corner. Your new report will appear under Report History.

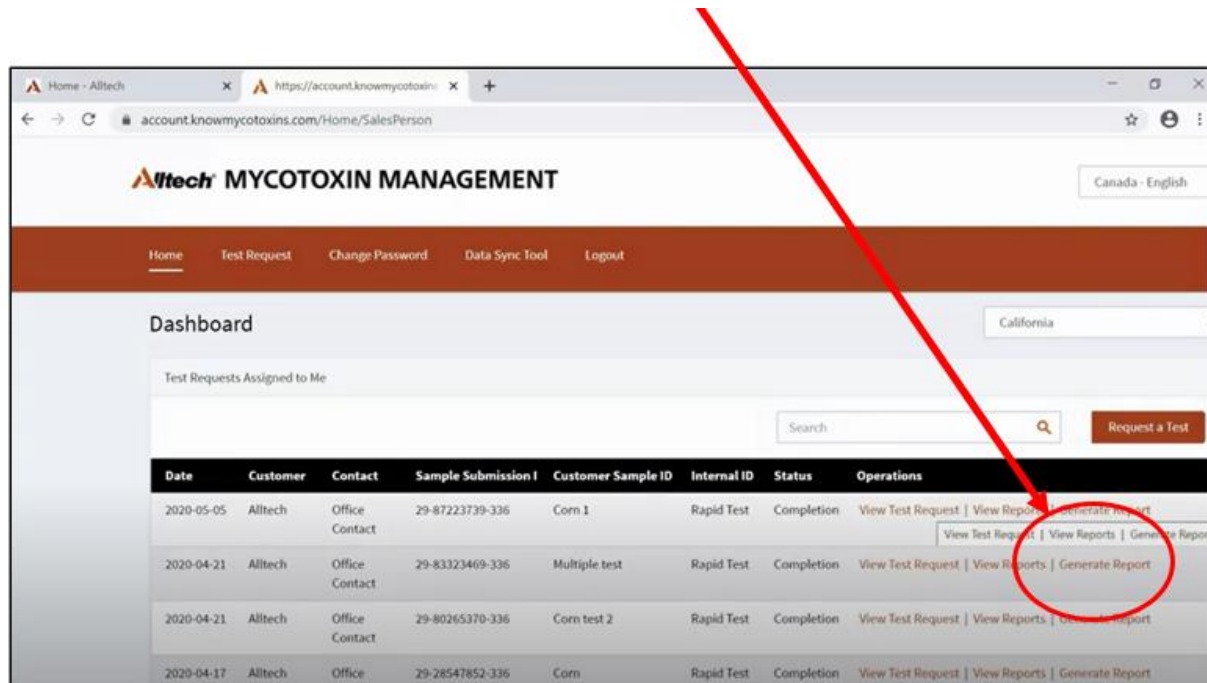


Step 5. Viewing a mycotoxin quick report on the Alltech RAPIREAD app


- A. Open the Alltech RAPIREAD App and login if required.
- B. Click on the **Submission** tab at the bottom of the screen.

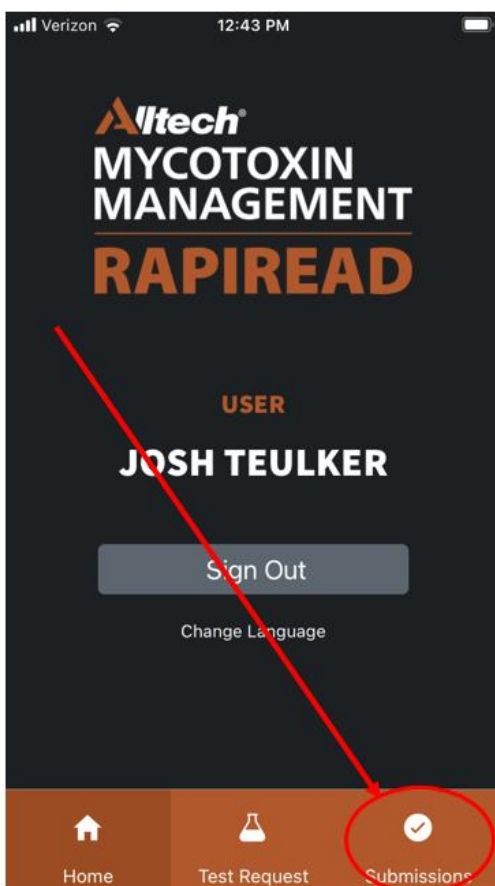


C. From here, you can view a full list of all your mycotoxin test submissions.

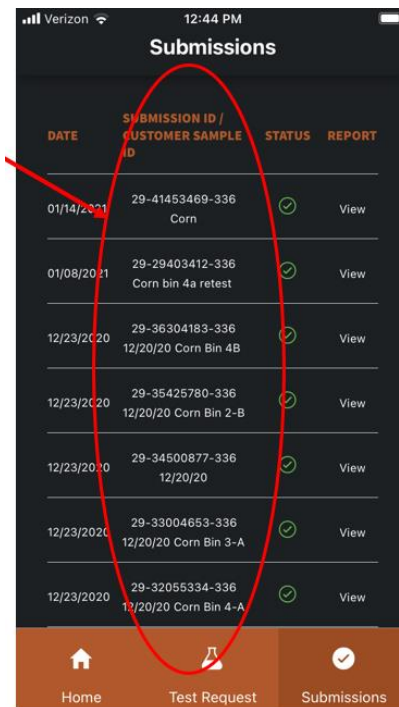

D. Under the report column, click **View** to view the results of a mycotoxin test.

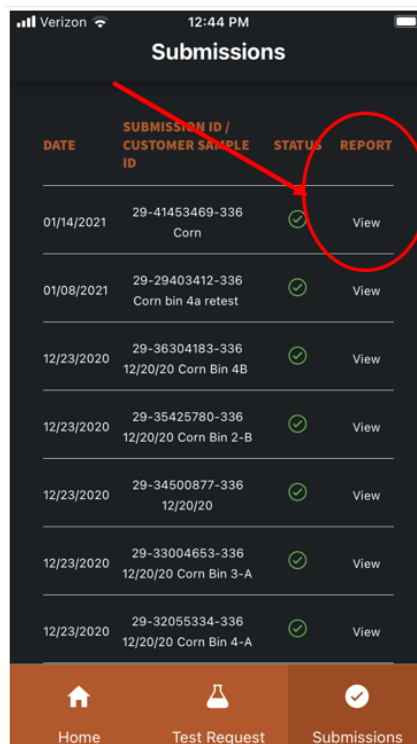

- E. Note: This can only be done when the submission from the App and the results from the Raptor have been synced. If the status column reads 'pending', you will need to sync the App and results.
- F. Click **Back** to navigate back to the view of all your test submissions.

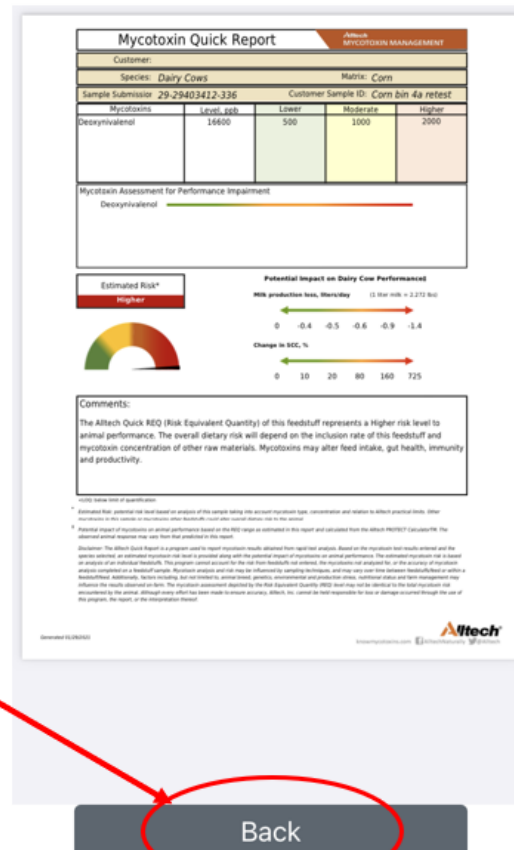

Overview video and further questions

You can view a short video at this [link](#) which outlines the main steps required to submit a mycotoxin test, run the test on the Raptor device and view the result on your portal and Alltech RAPIREAD app.

If you have further questions on getting set up or operating Alltech RAPIREAD please contact Josh Teulker at jteulker@alltech.com.